

TOTAL SULFUR AND TOTAL NITROGEN ANALYZER HELIOS S&N

ACCORDING STANDARDS

ASTM D5453 / ASTM D6667
ASTM D7183 / ASTM D7551
UOP 987 / IP490 / SH/T 0689
ISO 20 846

ASTM D4629 / ASTM D6069
ASTM D5762 / ASTM D7184
UOP 981 / UOP 971 / SH/T 0657
ISO 11905



 **TotalEnergies LICENSE**


s Associés
The Elemental Analysis



contact@eraly.com

HELIOS S&N

HELIOS S&N is the most sensitive and accurate analyzer for sulfur and/or nitrogen measurement dedicated to Refineries, Quality Control as well as contract laboratories.

It is used to measure rapidly Sulfur and/or nitrogen in liquid, oil, biofuel, hydrocarbon, solvent, lubricant oil, water, gas, LPG and solid samples.

Developed in collaboration between ERALY and end users (laboratory technicians and researchers) this equipment offers unrivaled advantages in ergonomics, reliability and robustness.

HELIOS S&N analyzer is the combination of 2 detectors (Sulfur & Nitrogen), without any alteration of their individual performances.

IT'S IMPORTANT TO HAVE CHOICE

For liquid samples:

- AutoInjector 1 position (single shot).
- Mini AutoSampler 6 positions, and rinsing vial plus trash.
- AutoSampler 50 positions, with several options (cooling, heating...).

For gas and LPG samples:

- AutoSampler for both gases and LPG (even mixtures).
- AutoSampler only for LPG (even mixtures).
- AutoSampler only for gases.

For calibrations:

Thanks to the linear response of the detector, it is possible to make,

- Calibration multipoint with 3 certified samples or more (calibration curve).
- Calibration one point (with single certified sample).

For configurations:

- Analyzer with separated computer.
- Analyzer with touch screen.
- Horizontal furnace.
- Vertical furnace.

Our analyzer HELIOS S&N is the most stable, it has a very long term stability with no deviation of the calibration during several months.

The Software HELIO-SOFT export function allows to calibrate with single one or more certified samples, in solid, liquid, gas or LPG.

No matter the type of matrix used for calibration, it is possible to convert the coefficient and use it to measure sulfur and/or nitrogen in all types of matrix.

It is possible to make various calibrations, according measurement ranges as well as method of calibration and save them in the software.

NEW GENERATION DETECTORS - Sulfur and Nitrogen



Thanks to the configuration Back-end sensor this new generation of **Sulfur detector** is more sensitive and **up to 20% more accurate** than the previous.

The **Nitrogen detector** has an on/off function of measuring compensation of chamber pressure, sample pressure, internal temperature in accordance with the values during calibration. This function among others, helps to increase accuracy of measurements.

Ultra low power consumption:

an environmentally-friendly and cost-saving analyzer. Pressure and temperature compensation for high measurement accuracy. Breakthrough design for power saving as well as thermal insulation.

Animated synoptic, with Touch screen 7":

at a glance, see, intensity of the UV Lamp signal, intensity of the PM signal; pressure of the measurement chamber, internal temperature, lifetime of filters.....

6 languages:

French, English, Spanish, German, Chinese, Russian.

Sulfur Detector



Nitrogen Detector



Service assistance inside:

detects early signs of trouble, allows predictive maintenance, identifies the needed service. All alarms could be displayed directly on the detector screen. On this screen you can display a picture of the concerned component and/or area.

For each trouble a table will show the minimum and the maximum accepted values, as well as the real time value.

Thanks to the online checking and assistance we can handle remotely the entire device.

WHY CHOOSE HELIOS S&N



- Acquiring two analyzers in single frame, at a price comparable to that of a single analyzer.
- Reduced costs for consumables. Commons parts: furnace, gases, electrical interfaces.
- Save time during your measurements, dosing of one or two elements in one injection.
- Reduction of operator manipulations and save space in your laboratory.
- Reduced maintenance costs, reduced maintenance time.

TECHNICAL SPECIFICATIONS

Injections	20 to 80 µl liquids / 10 ml gas / 20µL LPG		
Supply gas	Oxygen 99.998% - 3 bar	Argon 99.995 % - 3 bar.	
Electric	110 V / 230 V / 60Hz - 50 Hz / 800W		
Furnace	Temperature Furnace F2 up to 1100°C		
Dimensions (W x H x D)	51 cm x 83 cm x 56 cm / weight 32 kg		
SULFUR DETECTOR			
Detection Limit / Quantification Limit	5 ppb to 10%	/	10 ppb to 10%.
Accuracy RSD %	At 0.1 ppm 0.48 %	At 1 ppm 0.32 %	At 100 ppm 0.84 %.
NITROGEN DETECTOR			
Detection Limit / Quantification Limit	5 ppb to 10%	/	10 ppb to 10%.
Accuracy RSD %	At 0.1 ppm 0.51 %	At 1 ppm 0.40 %	At 100 ppm 0.88 %.

Analyzer with vertical furnace configuration and analyzer with horizontal configuration furnace has the same performances mainly for liquid and gas samples.

PYRO-COMBUSTION SYSTEM




As a manufacturer of tubular furnaces for industry and scientific research, we pay particular attention to the design and manufacturing of the combustion systems and combustion tube for our analyzers.

Nowadays, some famous international brands use our furnaces for their analyzers, this is the sign and proof that we have wide know-how and better technology.

The combustion efficiency, has a direct impact to measurement results. Combustion with injection and oxydation temperature well controlled, without soot production is one of the key of reliable measurement.

- Very robust furnace, guaranteed up to 5 years with no extra cost.
- Designed to facilitate quick replacement of thermocouple and combustion tube.
- Very low energy consumption with temperature accuracy better than +/- 1°C.
- No need to use gases (inert and oxygen) during heating, before reaching working temperature.

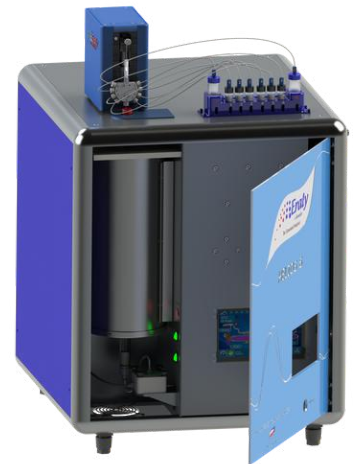


NEW SOFTWARE HELIO-SOFT



User-friendly and intuitive software.

Function export results to LIMS, to Excel and others commonly used formats such as PDF, HTML, XLS, CSV, or TXT.

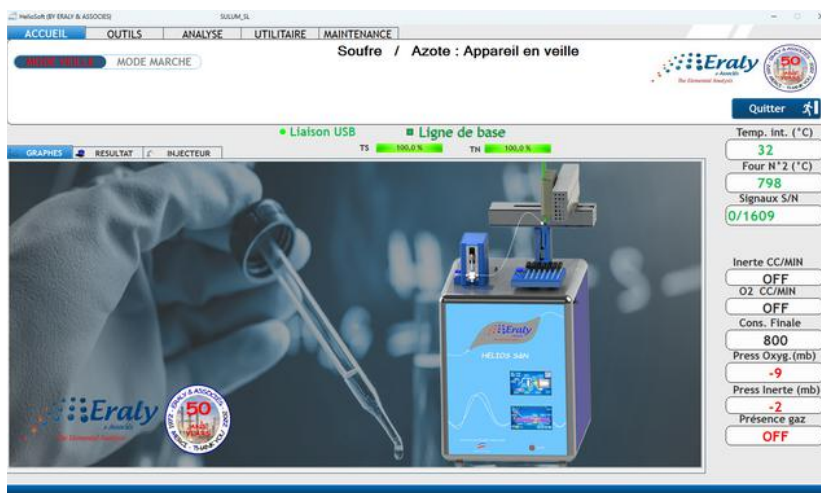

- Verification of calibration stability with graph (type control card).
- Alert function for tracking and replacement of spare parts and consumables.


ACCESSORIES



Mini AutoSampler / Autoljector

Tray of 1 position (single shot) or 6 positions, plus trash can and rinsing bottle.

Used for sampling and injecting liquids samples (e.g: methanol, petrolether, ethanol, toluene, naphtha, additives, kerosene, gasoline, diesel, bio diesel, bio ethanol, lubricant oil, waxes, organic solvents.....)

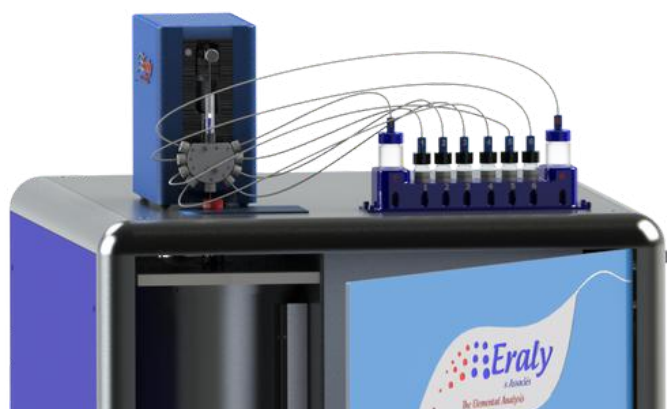

Gas and LPG SAMPLER

Compatible with standard: ASTM D6667.

Composed of 2 sampling loops, one for gas and one for LPG.

Used for sampling and injecting Butane, Propane, Methane, Natural Gas, Hydrocarbons and their mixtures (ethylene, methane, calibration gas mixes, etc.).

AutoSampler - 50 Positions

Tray of 50 positions, with cooling and/or heating options.

Used for sampling and injecting liquids samples (e.g: methanol, petrolether, ethanol, toluene, naphtha, additives, kerosene, gasoline, diesel, bio oil, lubricant oil, waxes, organic solvents.....).

Sampling volume adjustable between 10 μ l and 80 μ l. Sampling speed adjustable.

